152351

**Zoom-Stereomikroskop SVZB / SVZT**

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**Stereo Zoom Microscope SVZB / SVZT**

Operating instructions 4

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### Operating instructions

#### Introduction

In purchasing a NERIOX Stereo Zoom Microscope, you have chosen a quality product. It is suitable for use in schools and laboratories as well as in industry.

The Stereo Zoom Microscope consists of two separate microscope tubes, which are combined into one unit to allow simultaneous focusing on the object. Each tube has a prism, an achromatic lens and a wide field eyepiece. This creates a large, flat field of view. Both eyes view the object at a slightly different angle, which creates a stereoscopic effect.

With normal use, the maintenance requirement is reduced to a minimum.

These operating instructions provide information on the design, handling and maintenance of the microscope.

#### Design of the Stereo Zoom Microscope

Below you will find the designation of the individual components, which are also shown in the illustration:

1. Eyepiece
2. Dioptre correction
3. Prism housing
4. Zoom adjustment (2 rotary knobs)
5. Lens housing
6. Incident illumination
7. Rotary knob transmitted light
8. Rotary knob incident light
9. Main switch
10. Object clamp
11. Object plate
12. Stand base with transmitted light illumination
13. Head support
14. Locking screw (3 pcs.)
15. Focusing (2 rotary knobs)

#### Operating principle of the individual components

The Stereo Zoom Microscope consists of a stand with a head support, into which the stereo head is inserted and then fixed with the locking screws. Always hold the microscope by the stand when you want to move it.

#### Tube

The head of the Stereo Zoom Microscope is equipped with a 45° tube that can be rotated through 360°. The head also has dioptre correction on both eyepiece tubes.

#### Optical specifications (versions)

The Stereo Zoom Microscope has 2 zoom lenses and a pair of wide field WF10x eyepieces.

Model no.	Eyepieces	Lenses	Magnification	Version	Tube
152351.0100	WF10x	0.7x ... 4.5x	7x ... 45x	binocular	45°
152351.0200	WF10x	0.7x ... 4.5x	7x ... 45x	trinocular	45°

#### Object stage

The stand is equipped with an object stage with a transparent object plate and two object clamps. A black and white object plate is provided as a standard accessory.

#### Focusing

Focusing is used to bring the image of the object into focus.

### Adjusting the magnification

The Stereo Zoom Microscope has 2 achromatic zoom lenses. The magnification is continuously adjusted by turning the rotary knobs for the zoom function.

### Illumination

The stand has adjustable incident light and transmitted light LED illumination. The main switch is located on the right side of the stand base. The two lights can be used simultaneously. The rotary controls for incident light and transmitted light are located on the right side of the stand base.

- LED: Incident light illumination
- LED: Transmitted light illumination

### Preparing the Stereo Zoom Microscope for use

Remove the microscope from its packaging and place it on a stable table. Insert the mains plug into a power socket and operate the main switch. Then switch on the illumination of the microscope. Then slide the eyecups over the eyepieces. Now sit comfortably in front of the microscope and look through the eyepieces.

### Attention:

When using the black and white object plate, the transmitted light illumination must be switched off!

### Working with the Stereo Microscope

- Set the focus to a neutral position.
- Place an object on the object plate and adjust the height of the stereo head until you see a reasonably sharp image and fix the head support with the locking button on the rear side.
- Move the dioptre correction from both eyepieces into the neutral position (marking groove on the tube).
- Adjust the pupil distance by moving the prism housings towards or away from each other until you see a round image with both eyes.

### Adjusting the eyepieces to get a sharp image over the entire zoom range

- Set the magnification to the largest value using the zoom adjustment function. Now focus on a specimen that is as flat as possible (e.g. a text).
- Set the magnification to the smallest value.
- Now correct the sharpness of the left eye by closing the right eye and adjusting the dioptre correction. Do the same with the right eye.

### Maintenance and cleaning

Always cover the microscope with a dust cover after use. Keep the eyepieces mounted in the microscope tube at all times, to prevent dust from getting into the device.

### Cleaning the optics

If the lenses of the eyepiece are dirty, they can be wiped with a piece of lens paper (circular movements). If this does not bring an improvement, moisten the lens paper with a drop of xylene or alcohol. Never apply xylene or alcohol directly on the lenses!

If dirt is clearly visible in the field of view, this is located on the lowest lens of the eyepiece. The eyepiece can simply be pulled out of the tube. Now clean the outside of the lens. If dust is still visible, make certain that the dust is really in the eyepiece by rotating the eyepiece. If this is definitely the case, carefully unscrew the bottom lens from the housing and clean it.

### Warning:

Cloths with plastic fibres can damage the coating of the lenses!

### Cleaning the stand

Dust can be removed with a brush. If the stand is very dirty, the surface can be cleaned with a non-aggressive cleaning agent.

### Procedure for illumination problems (incident light / transmitted light)

The built-in LEDs have a very long service life. If there is a problem with the illumination, please send the microscope to our Service Center → [Servicecenter@brw.ch](mailto:Servicecenter@brw.ch)

### Warning:

LEDs or electronic parts may only be repaired or replaced by qualified personnel! In the event of non-compliance, any liability will be refused.

### Optional accessories

152371.0010	Eyepiece (pair)	WF15x / 15
152371.0020	Eyepiece (pair)	WF20x / 10
152371.0030	Eyepiece (pair)	with reticle (1 piece) WF10x
152371.0050	Attachment lens	2.0x

### Repairs

If your microscope needs to be repaired, please contact our Service Center → [Servicecenter@brw.ch](mailto:Servicecenter@brw.ch)

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